




# ***SureQuant: Novel Turnkey Targeted Toolbox for Routine Quantitative Pathway Proteomics***

Aaron Gajadhar, PhD  
Strategic Marketing Scientist – Quantitative Proteomics


# Fundamental Challenge of Biology is to Reliably Quantify Protein Components

## Classical, mechanism-focused protein analysis

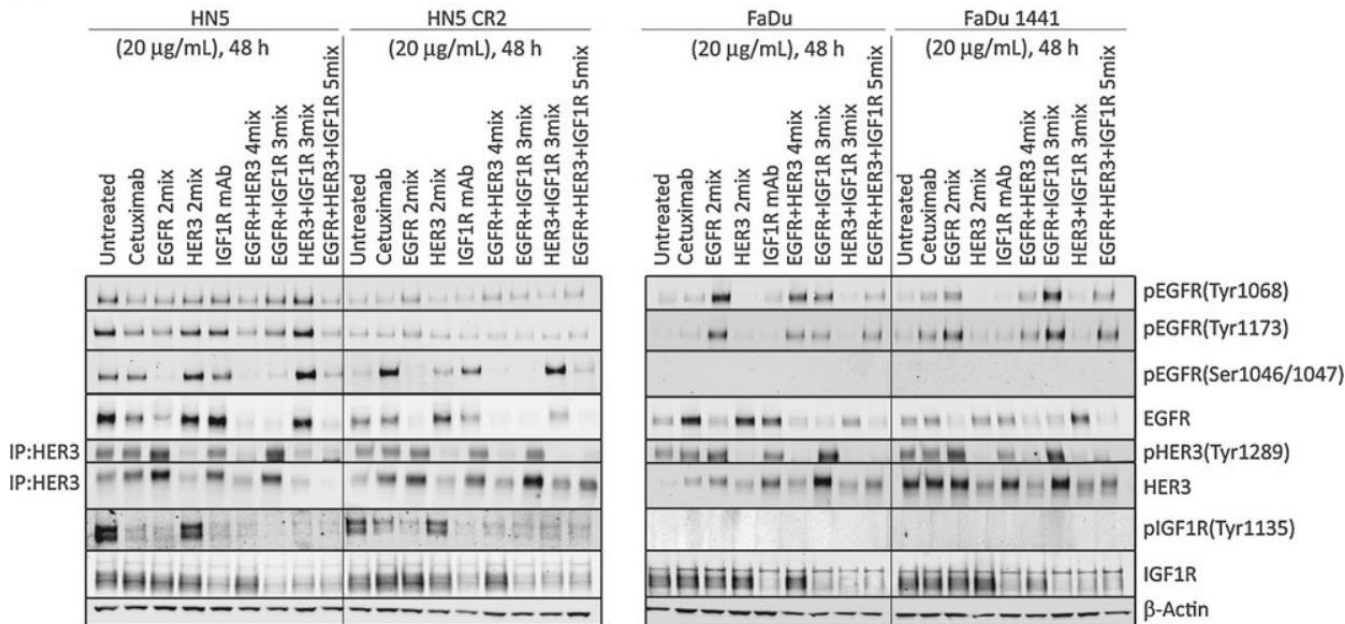

*Mol Cancer Ther.* 2016 Jul;15(7)

## Systematic proteomic characterization

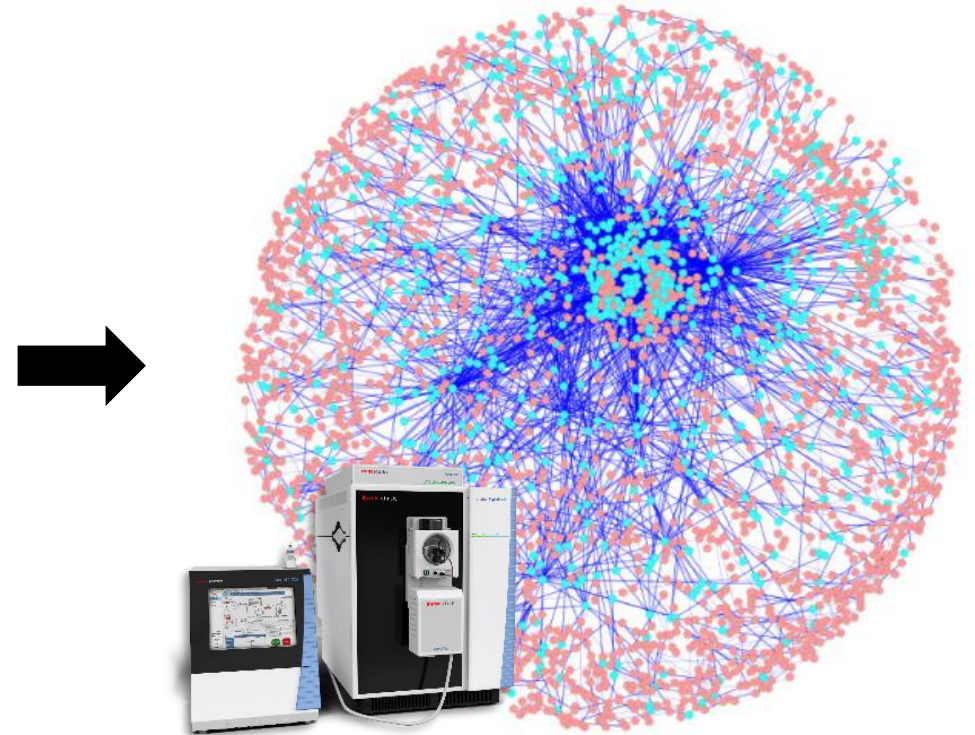

*BMC Syst Biol.* 2011 Aug 5;5:122

**Need to know how the proteome profile is changing: Time, disease, stimulus, environment, ...**



## Tymora Developing Extracellular Vesicle-Based Phosphoproteomic Liquid Biopsies

## Early Results From Second Side-Out Study Indicate Benefits of Molecularly Guided Cancer Treatment

## MSK Researchers Develop High-Sensitivity Mass Spec Assay for Clinical Cancer Work

## Molecular Cell

## A Strategy to Combine Sample Multiplexing with Targeted Proteomics Assays for High-Throughput Protein Signature Characterization

Technology

“...profile **tumor signaling pathways** via blood or urine samples, similar to how liquid biopsies make use of ctDNA analysis...”

“...number of research groups are exploring the use of **phosphoproteomics** analysis to aid selection of patient treatments and track their response to therapy.”

“...using **molecular data to guide patient treatment** extends progression-free survival in metastatic breast cancer patients.”

“...bolster the case for **phosphoproteomic** data specifically, as this was the primary type of data used for guiding therapy.”

“ We are working on clinical tests to measure **signaling in cancer specimens** as part of diagnostic clinical tests...”

“...more **quantitative measurements than antibody assays** and the ability to **profile pathways more broadly**, covering molecules for which there are not good antibodies.”

“ We **targeted 131 peptides** corresponding to 69 **cancer-related proteins**

“ We utilized TOMAHAQ to assay expression of 69 protein targets across the **entire NCI-60**

“ BAZ1B, also a tyrosine-kinase....expression was shown to correlate with **drug response sensitivity**...”

# Quantitative LC-MS for Pathway Proteomics - Promises & Pitfalls



**Extreme sensitivity and inherent specificity**



**Quantitate 100s - 1000s of proteins**



**Identify PTMs and protein interactions**




**Compatible with enrichment and isolation methods**



**Detect protein isoforms and modifications**



## LC-MS Pathway Proteomics

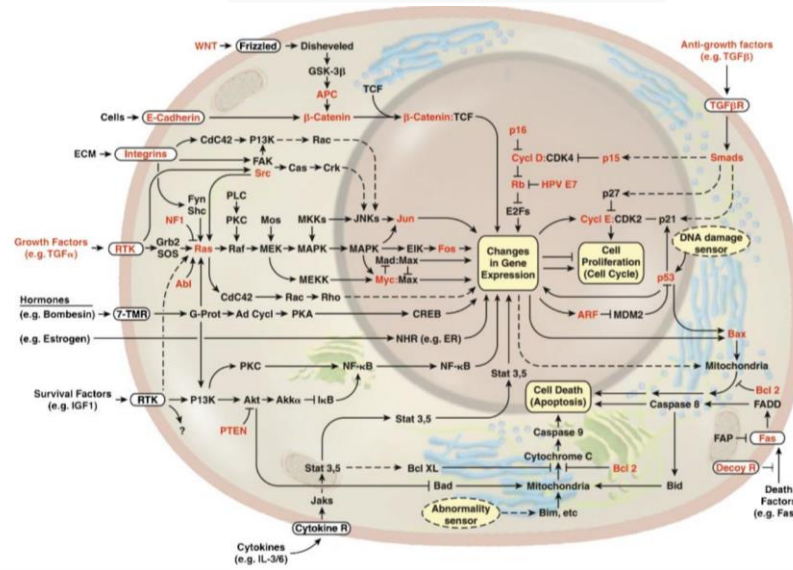

Cell. 2000 Jan 7;100



**Dynamic range of target abundance**



**Quantitative reproducibility 'Missing values'**



**Tradeoffs for discovery vs. targeted workflows**



**Tedious, time-consuming assay development**



**Lack of truly integrated and standardized workflows**



# SureQuant Targeted Mass Spec Assay Kits


Validated, modular reagents for multiplexed target protein quantitation



Thermo Scientific™ SureQuant™ Mass Spec Assay Kits

## Highlights

- **Complete**—includes all reagents for successful monitoring of MS system performance, sample preparation, and absolute quantitation of target proteins/peptides
- **Verified**—antibodies, peptides, and control lysate are rigorously tested for specificity and successful quantitation of each target peptide
- **Multiplex**—able to quantitate different targets (including isoforms and phosphorylated states) from the AKT/mTOR pathway
- **Flexible**—modular format allows for immune-enrichment only, or in combination with relative or absolute quantitation

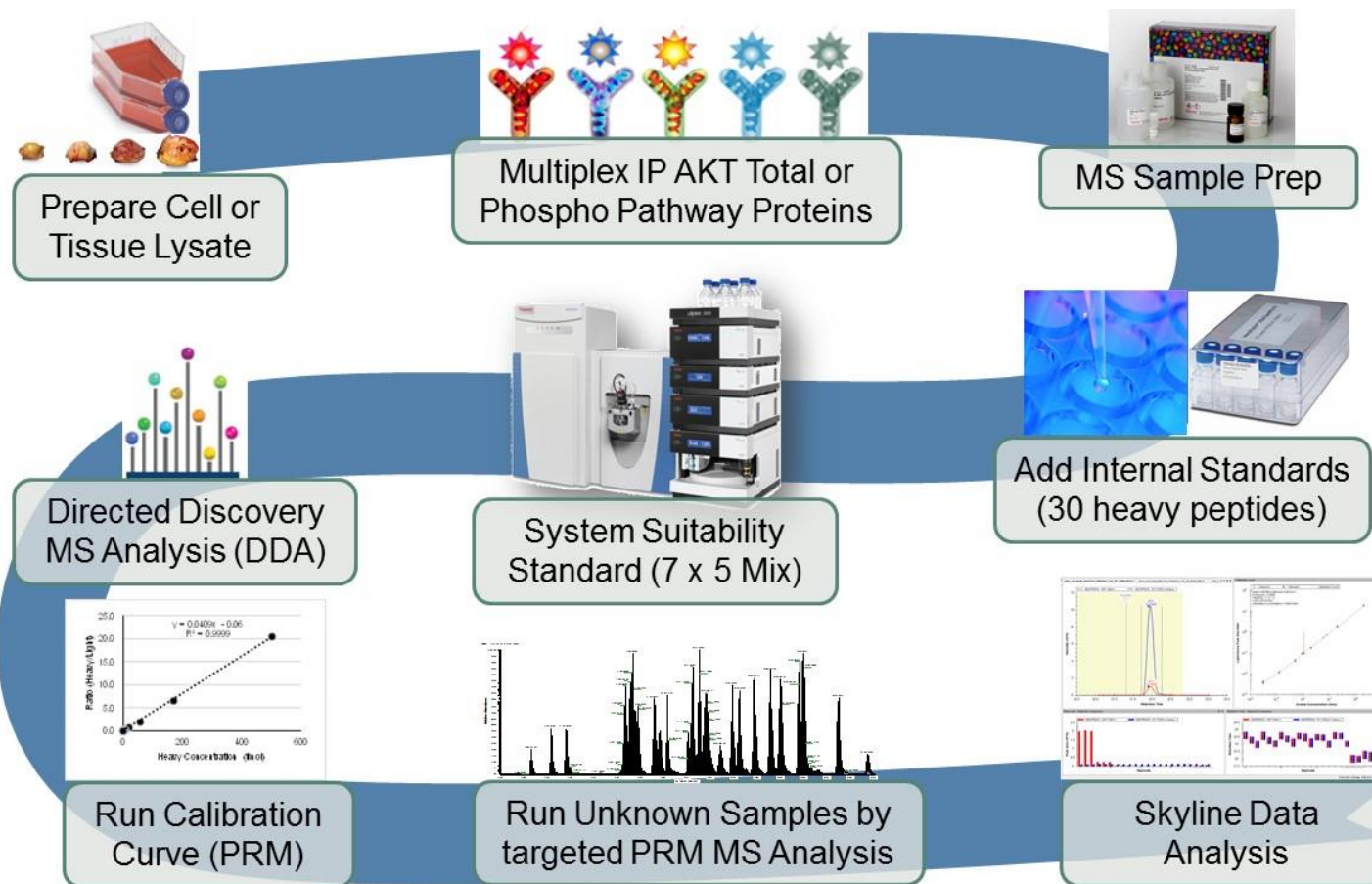

# The Next Leap Towards Turnkey Pathway Quantitation

Two new 'data aware' intelligence-driven systems to meet the evolving needs of the protein mass spectrometry community



**Thermo Scientific™ Orbitrap Exploris™ 480  
Mass Spectrometer**


*Robust workhorse for proteome  
analysis and quantitation*



**Thermo Scientific™ Orbitrap Eclipse™ Tribrid™  
Mass Spectrometer**

*Highest performance, flexibility, usability for  
protein structure and quantitation studies*


# Thermo Scientific™ SureQuant™ IS Targeted Protein Quantitation Solution

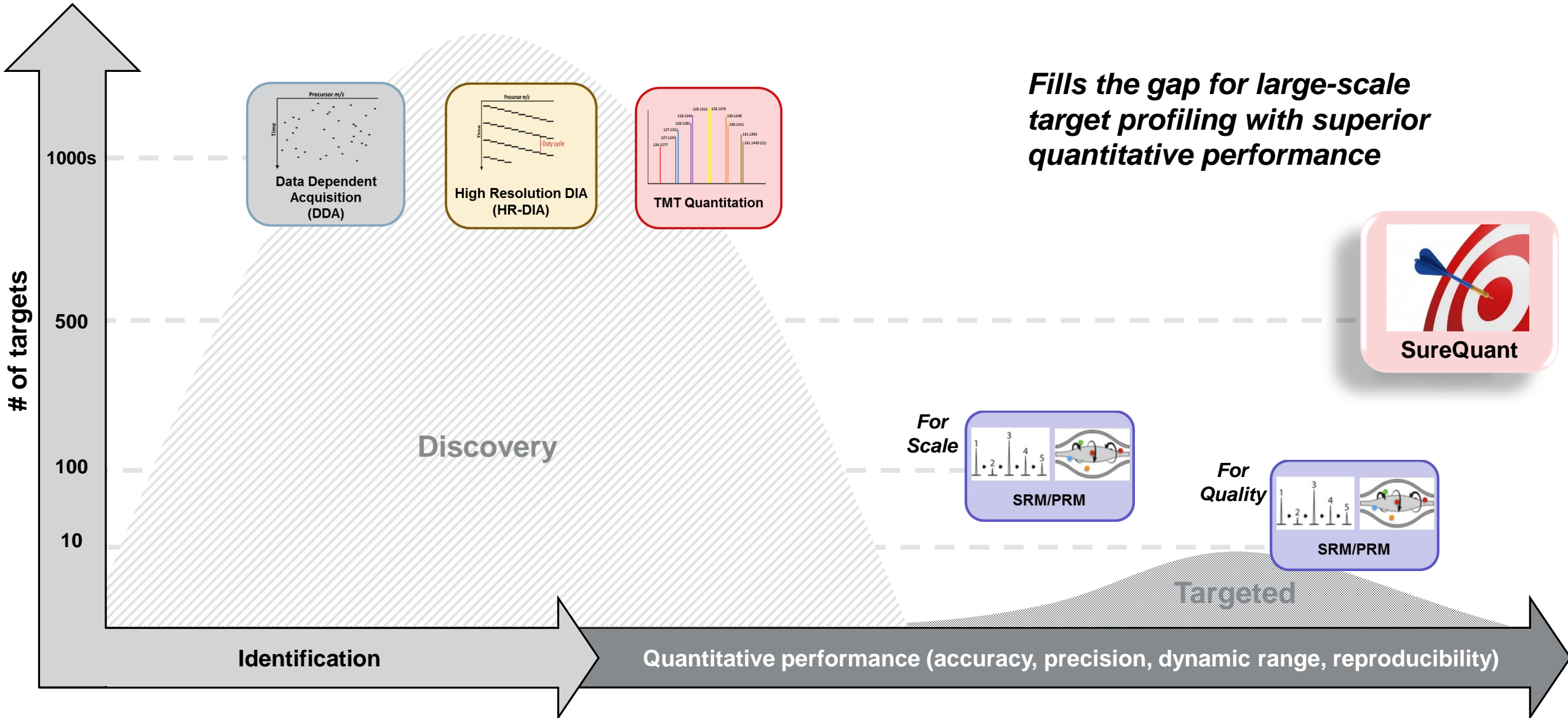

# Introducing SureQuant Targeted Protein Quantitation

Intelligent Data Acquisition Strategy for Sensitive, Precise and Reproducible Quantitation


Synthetic Heavy Peptides

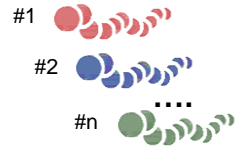



Sample



LC-MS Analysis

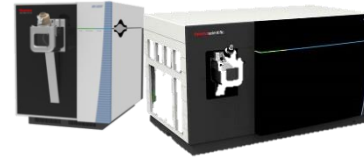

**SureQuant method**

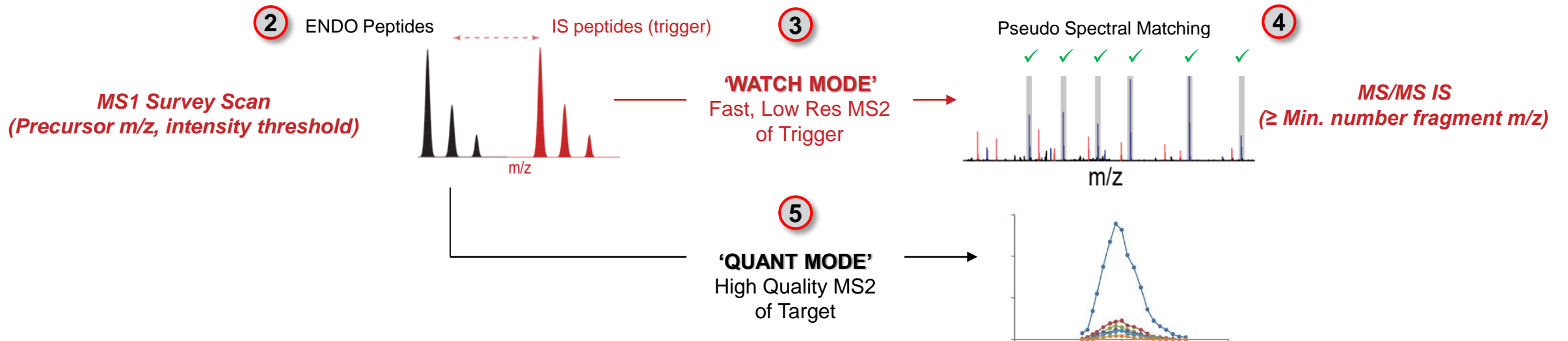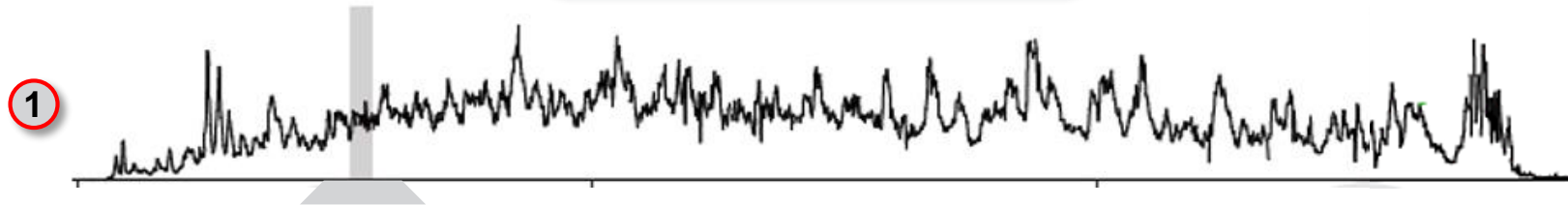



# Overall SureQuant Acquisition Workflow – Embedded Application Specific Kits

1

Survey Run

Ready-to-use template



LC-MS Analysis



Survey Run Analysis



Kit specific Skyline analysis template

Directed DDA

*Inclusion list optimal IS precursor ions*



Export reports



Targeted Mass

*Intensity response of precursor ion*

Targeted Mass Trigger

*'Optimal' fragments / selected peptide precursor*



2

SureQuant

Ready-to-use template

Internal standards



Sample



Import parameters



LC-MS Analysis



SureQuant



Target Quantification




Kit specific Skyline analysis template




# SureQuant Workflow – Validated Method Templates

Optimized Methods Enable Focus on Science not Set-Up

## System Template

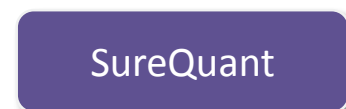





## Kit Specific method

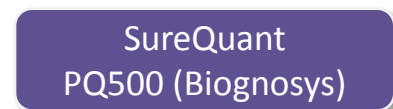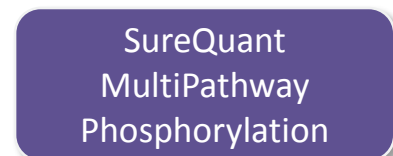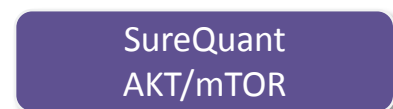

## Custom panel




Generic template

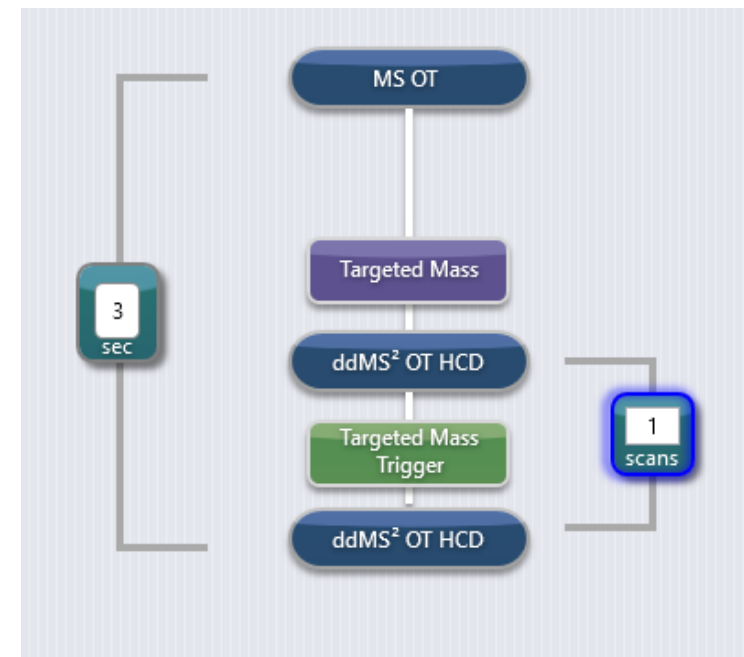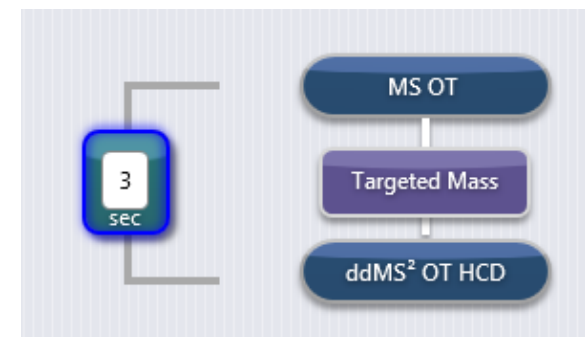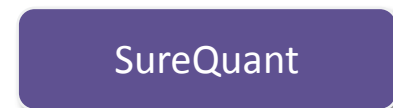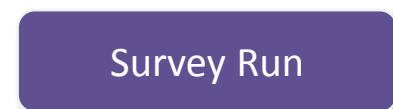

- Ready-to-use templates validated for application specific SureQuant peptide kits
- SureQuant available on all TNG Orbitrap instruments with Tune v3.3 or higher

# SureQuant Acquisition Drives Enhanced Productivity by Intelligent Detection of Targets


■ Non-productive MS2 scans    □ Productive MS2 scans

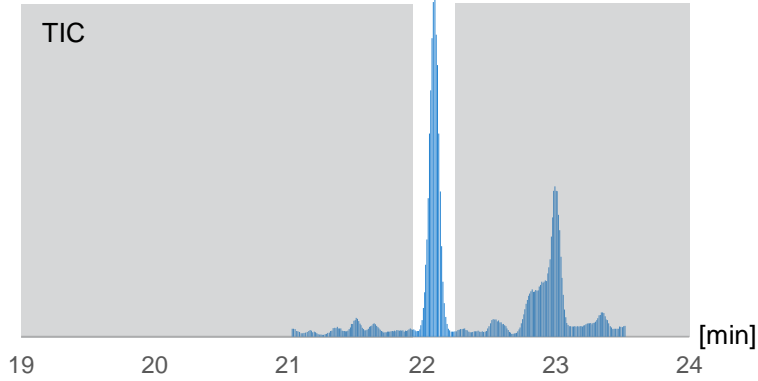


**PRM**

SIL


**LCDSGELVAIK**

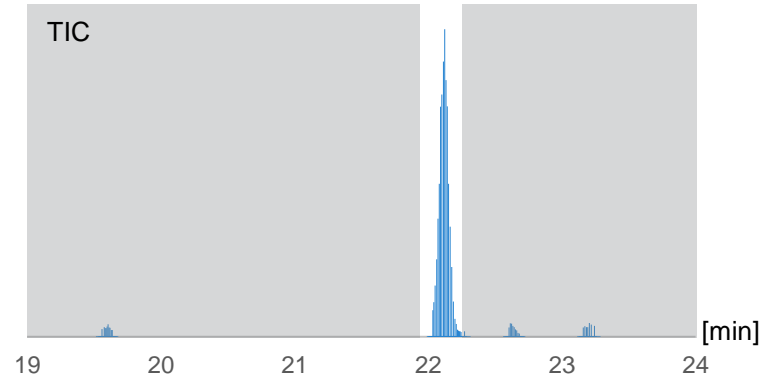

**SureQuant**

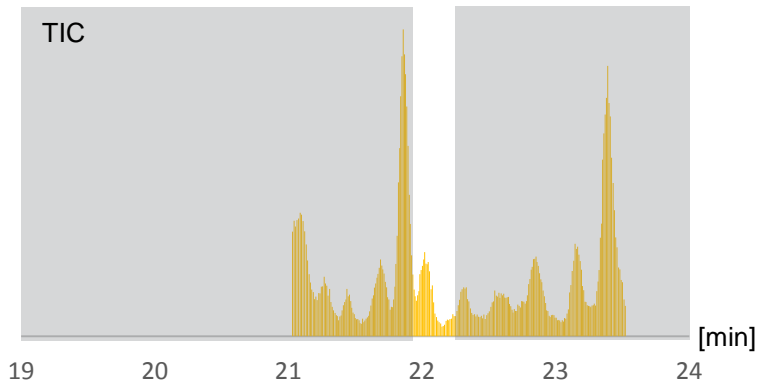

ENDO

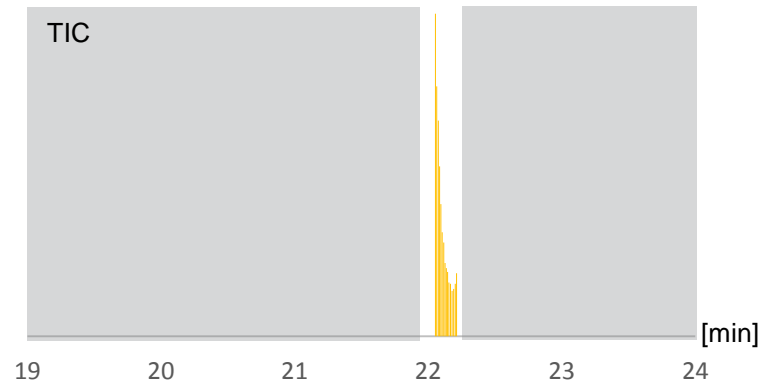

## Parameters / conditions

- PRM
  - 15k / 20 ms
  - 2.5-min monitoring window
- SureQuant
  - Watch mode: 7.5k / 10 ms
  - Quant mode: 60k / 116 ms
  - 10-ppm tolerance MS1 (w/ lock mass)
- Sample
  - 250 ng cell line digest
  - 50 fmol internal standards

**Proportion MS2 productive scans  
(% MS2 acquisition time)**

**PRM ≈ 10-15 %**




**SureQuant ≈ 80-90 %**

# SureQuant's Intelligent Acquisition Robustness Overcomes Analytical Fluctuations





- SureQuant target detection occurs independent of RT scheduling or poor chromatography conditions allowing reliable and robust target quantification

Original elution profile

5 min delayed elution profile

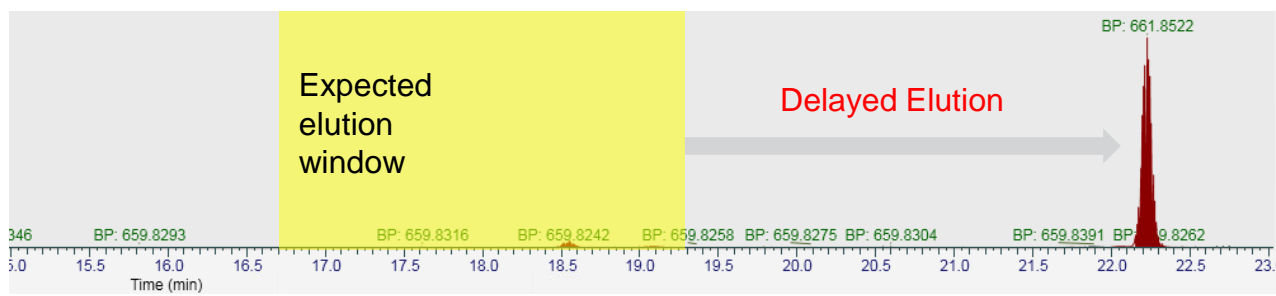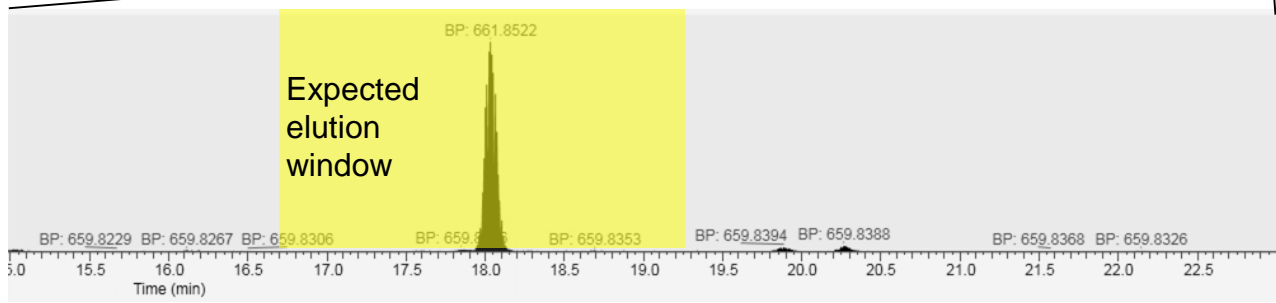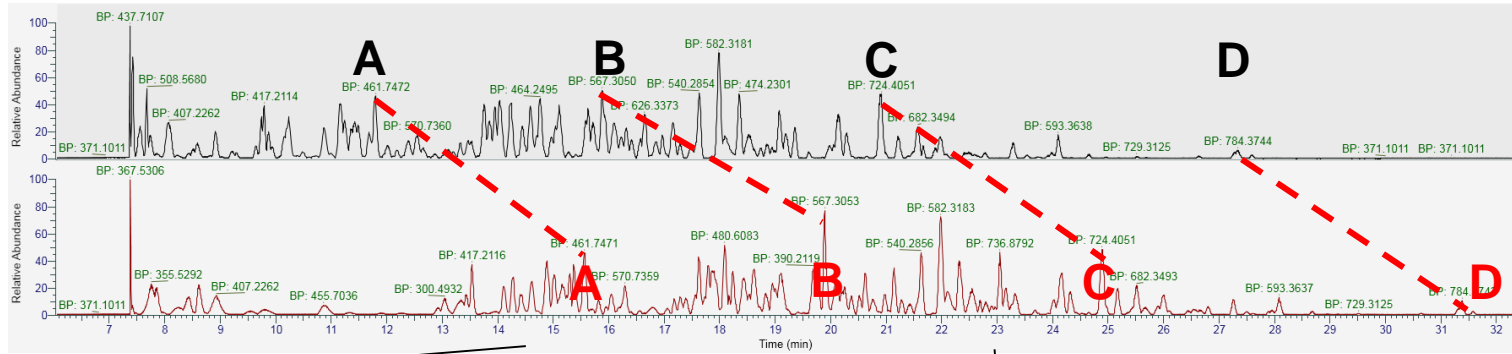

**Heavy Peptide Detection**

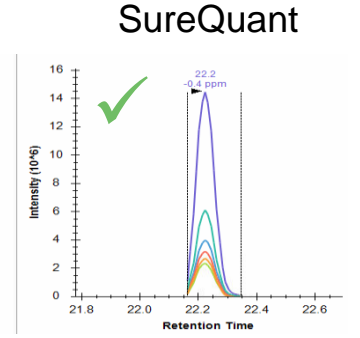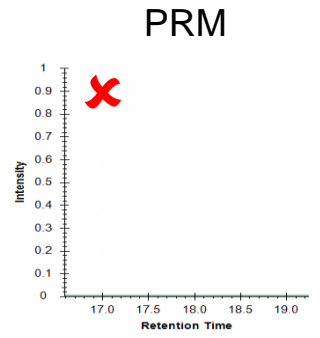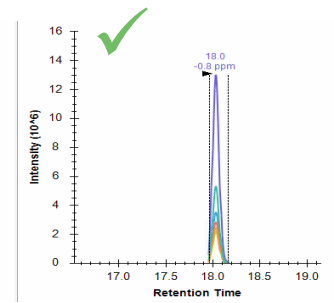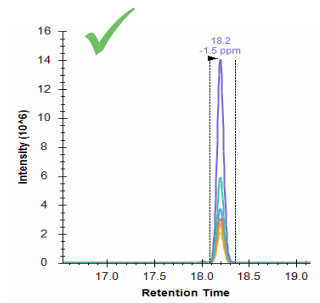




# Application Example #1 - AKT/mTOR pathway analysis

- Many genetic alterations in cancer cells modify the protein expression from AKT, RAS and TP53 pathways.
- Quantitative measurement of alterations in the expression of pathway proteins and post-translational modifications (PTM) is necessary for classifying disease states, monitoring cancer progression and determining treatment response.
- **SureQuant analysis of the AKT/mTOR pathway (12 protein/30 peptide targets)**
- **Capillary flow LC-MS setup (1.2 uL/min)**
- **Detection of targets without IP enrichment?**

## AKT/mTOR protein pathway

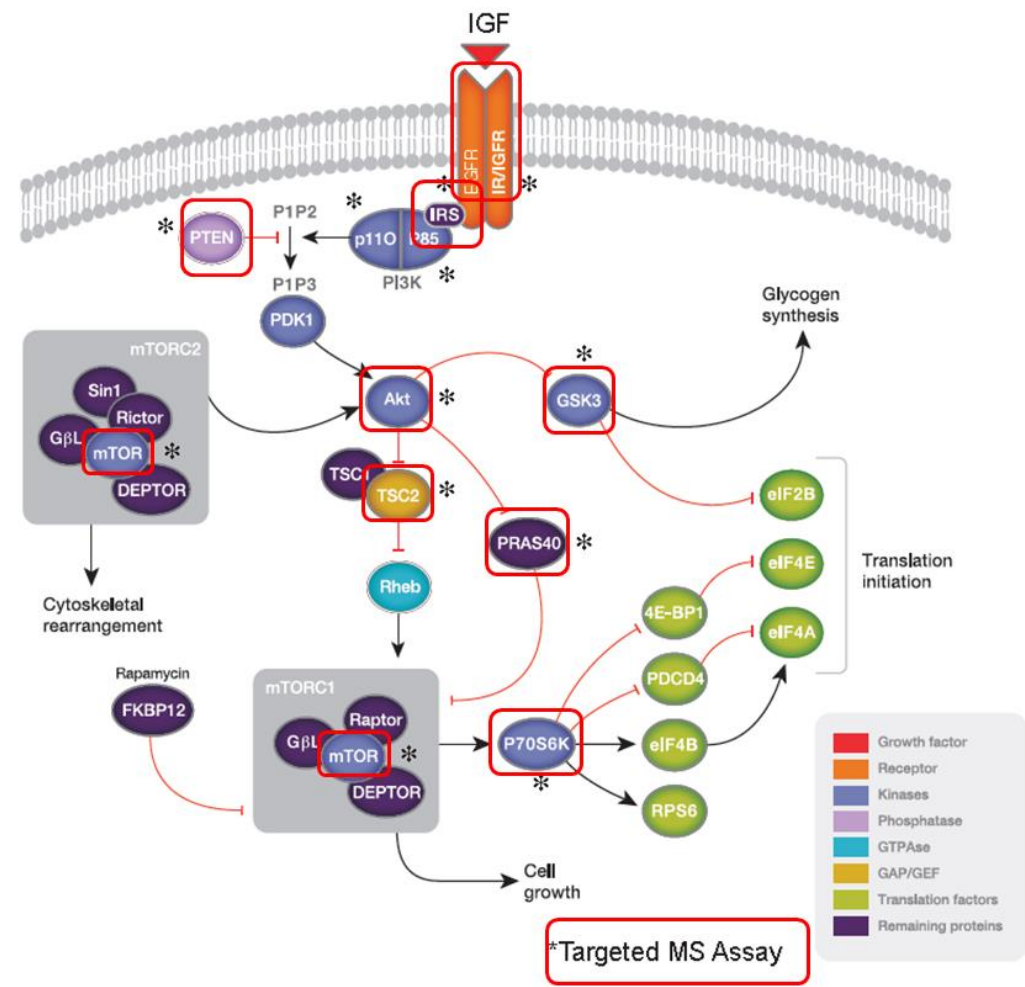

# SureQuant Allows Sensitive Detection of AKT/mTOR Proteins Without IP Enrichment

Protein	Peptide	PRM 15K-20ms	SureQuant 60K-116ms
GSK3β	TPPEAIALC[Carbamidomethyl]SR	1	20
AKT2	LPFYNQDHER	0	17
IGF1R	AENGGPGVGLVLR	1	15
PTEN	IYNLC[Carbamidomethyl]AER	0	15
GSK3β	LC[Carbamidomethyl]DSGELVAIK	5	15
GSK3β	LLEYTPRAR	0	15
GSK3α	VTTVVATLGQGPER	6	15
AKTS1	GALAEAAR	5	13
AKT1	NDGTFIGYK	0	13
AKT2	SDGSFIGYK	0	13
IGF1R	IDIHSC[Carbamidomethyl]NHEAEK	1	12
PTEN	AQEALDFYGEVR	10	11
MTOR	LFDAPEAPLPSR	11	11
MTOR	TLDQSPELR	2	10
AKT1	YSFQTHDR	0	10
KS6B1	FEISETSVNR	0	9
TSC2	SNPTDIYPSK	0	9
GSK3α	SQEVAYTDIK	1	9
TSC2	GQPEGPLPSSSPR	0	8
AKTS1	SSDEENGPPSPDLDR	0	8
IGF1R	TTINNEYNYR	0	7
MTOR	ETSFNQAYGR	2	6
MTOR	GNNLQDTLR	0	6
TSC2	GYTISDSAPSR	0	5
PTEN	NNIDDVVR	0	4
KS6B1	DGFYPAPDFR	0	1
AKT3	LVPPFKPQVTSETDTR	0	0
IRS1	SVSAPQQIINPIR	0	0
AKT3	TFHVDTPPEER	0	0
IRS1	TGIAAEEVSLPR	0	0

Number of peptides detected:


11

26

Number of spectra including ≥ 5 ref. fragments




30 min Gradient Method

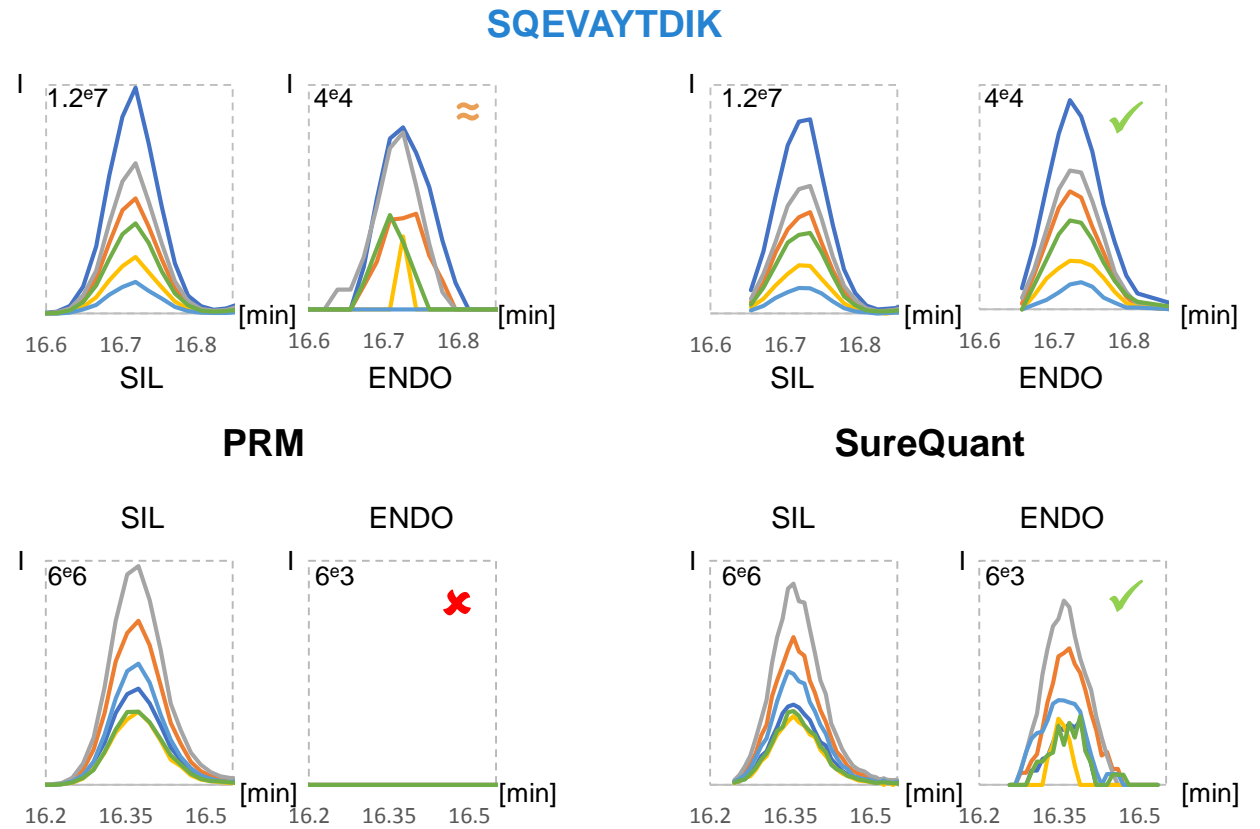

# Rapid Pathway Analysis – Sufficient Peak Sampling Ensures Precise Quantitation

- The entire pathway (12 proteins/30 peptides) was quantified with a **10 min gradient** (n=3 replicates)

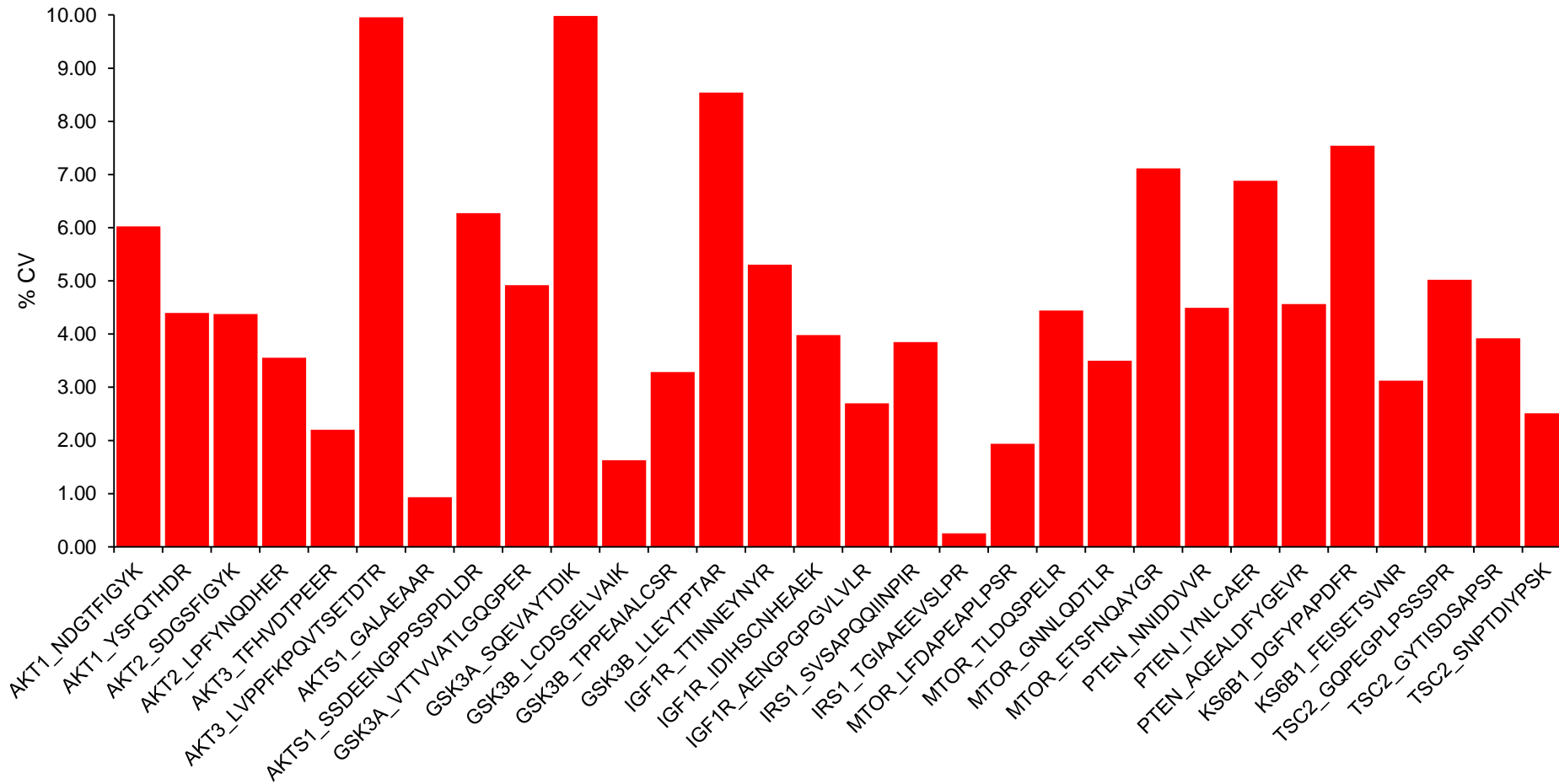

**100%**  
Peptides  
quantified  
<10% CV

**67%**  
Peptides  
quantified  
<5% CV

# Customer Application Example - Global pTyr Profiling In Human Tumor Specimens

Collaboration with  
Forest White Lab, MIT




*Tyrosine phosphorylation analysis  
of tumors uncovers druggable targets*


Tyrosine kinases  
**0.3%** of genome  
**30%** of oncoproteins



Human tumor

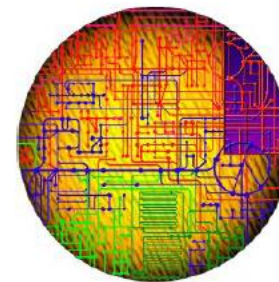

Signaling targets

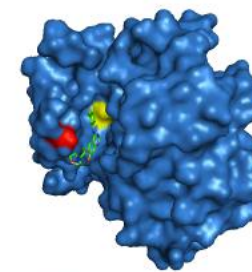

Therapeutics

## SureQuant triggered global pTyr profiling in colon tumor specimens

385 pTyr  
heavy  
peptide  
standards


Mix



Tumor specimen



pY peptide enrichment

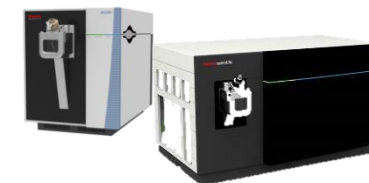


SureQuant Method

For research use only




# SureQuant pTyr Signaling Analysis Offers Reproducible and Complementary Insights


## Reproducible target quantitation

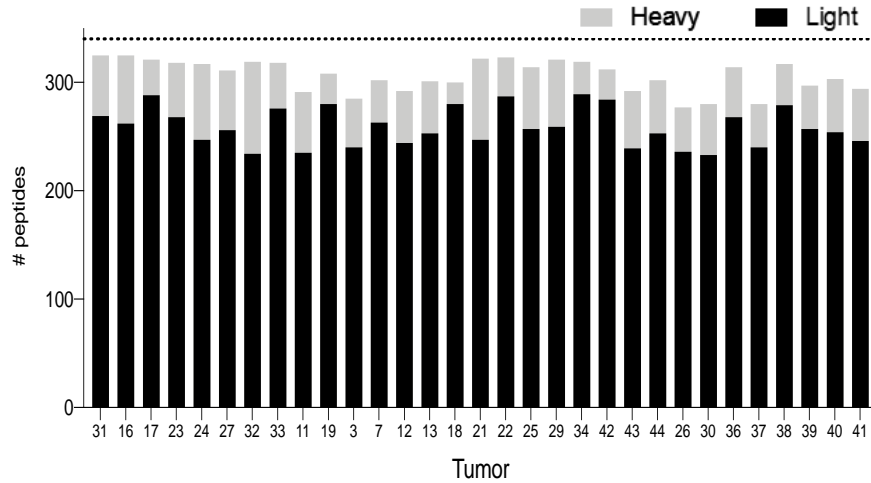

## Activation of actionable targets can be monitored across colon tumors

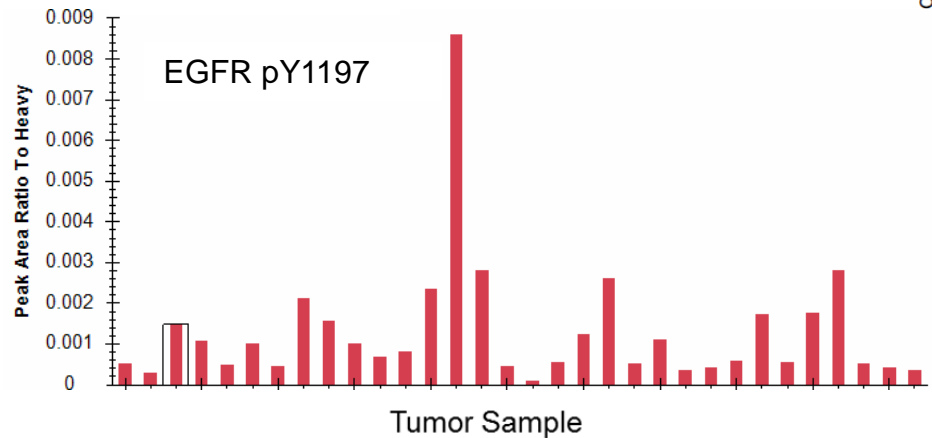

## Signaling heterogeneity uncovers biological insight not available from genetic subtyping

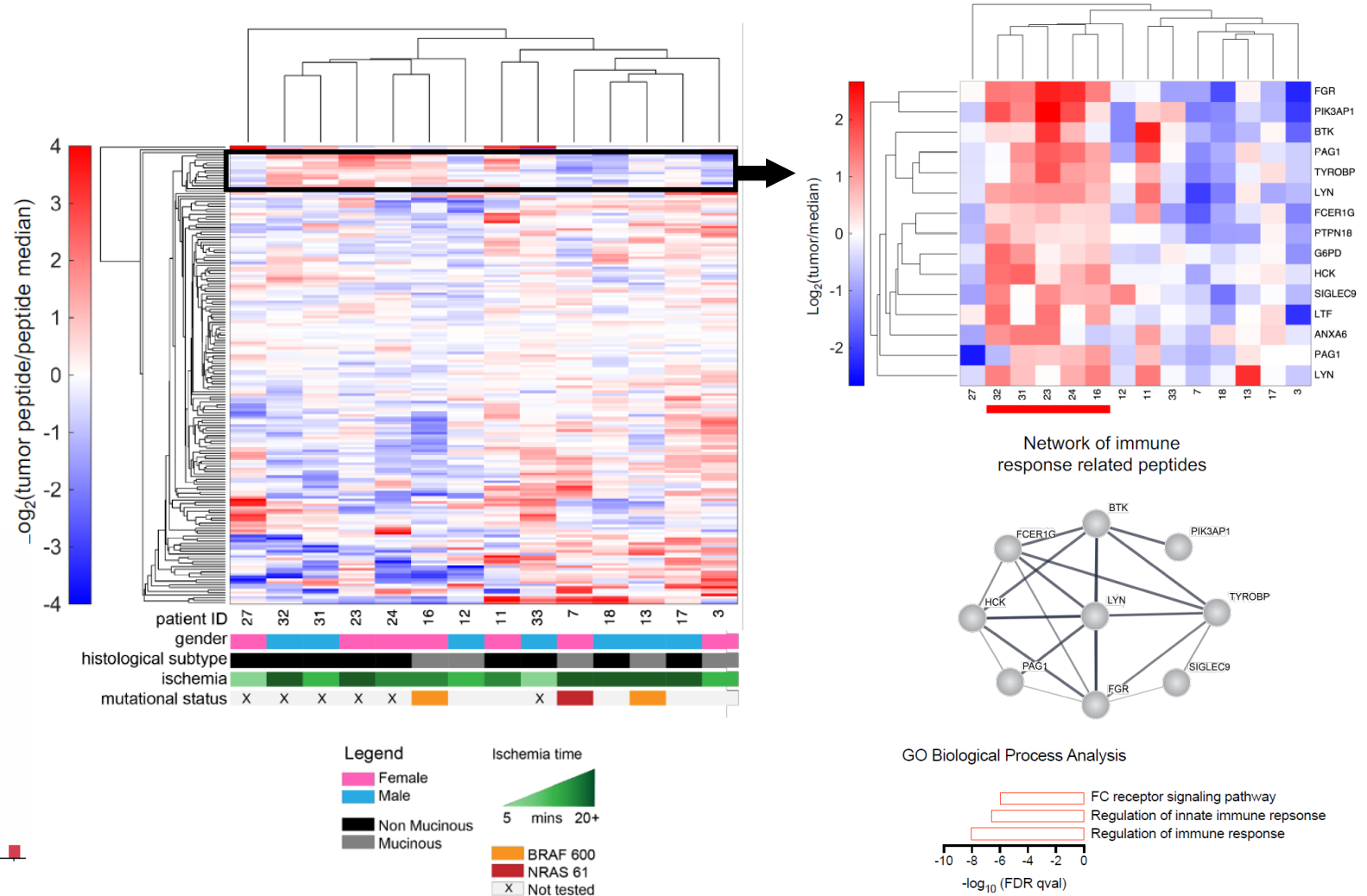

# SureQuant Large-Scale Target Profiling with Superior Quantitative Performance

SureQuant

PRM/SRM

PRM/SRM

up to 500 pairs of H/L targets/30 min

**Value Prop**

- \*Highest targeting efficiency
- \*Complete quan profile every analysis
- \*Maintain sensitivity *AND* target scale

\*Favoring sensitivity  
up to 70-80 pairs of H/L targets in 30 min  
(60k – 116ms)

\*Favoring target number  
up to 400-450 pairs of H/L targets in 30 min  
(7.5k – 10ms)

<b>Quantification Performance</b> (Precision/Accuracy)	★★★★	★★★★	★★★
<b>Sensitivity</b> (LLOQ)	★★★★	★★★★	★★★
<b>Scale</b> (# Targets)	★★★★	★★★	★★★★
<b>Efficiency</b> (Productive Scans)	★★★	★★★	★★★
<b>Load-and-play</b> (Minimal adjustment)	★★★	★★★	★★★

# Acknowledgements

70,000 Thermo Fisher employees in 50+ countries  
supporting the global business of science



## Applications Team

- **Sebastien Gallien**
- **Bhavin Patel**
- Khatereh Motamed
- Daniel Lopez-Ferrer
- Andreas Huhmer

## Product Management Team

- Tabiwang Arrey
- Christian Thoeing
- Kerstin Strupat
- Markus Kellmann
- Alexander Harder
- Derek Bailey
- Graeme McAlister
- Romain Huguet
- Shannon Eliuk
- Vlad Zabrouskov

## Marketing Team

- Julian Saba
- Sally Webb
- Maciej Bromirski
- Scott Peterman
- Lisa Thomas
- Ken Miller